Isothermal Master Mix for Genie®

製品コード: OP-IMM200, OP-IMM400

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OptiGene Limited
Unit 5 Blatchford Road Horsham
West Sussex RH13 5QR UK

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Isothermal Master Mix

For research purposes only. Not for use in diagnostic procedures for clinical purposes.

FOR *IN VITRO* USE ONLY

Product Description:

The Isothermal Master Mix is an optimised mastermix designed to simplify the preparation of an isothermal amplification assay. The mastermix contains a proprietary, fast, novel DNA polymerase, proprietary thermostable inorganic pyrophosphatase, optimised reaction buffer, Mg²⁺, dNTPs and a ds-DNA binding dye (FAM detection channel).

Stability:

Store the Isothermal Master Mix reagents at -20°C in a constant temperature freezer.

Usage:

A typical 25µl reaction would contain the following:

Component	Volume
Isothermal Master Mix	15.0 µl
Primers	5.0 µl
Template	5.0 µl
Total	25.0 µl

Typical amplification conditions for this kit are:

Amplification	30 mins	65°C
Anneal		98°C - 80°C
		- ramping at 0.05°C per second

Note:

Due to the presence of the inorganic pyrophosphatase there will be no accumulation of PPI.

We recommend the following Primer levels per 25µl reaction:

F3 & B3	5 pmol each primer
LoopF & LoopB	10 pmol each primer
FIP & BIP	20 pmol each primer

Quality Control:

The Isothermal Master Mix has been tested for performance in Loop-mediated Isothermal DNA amplification.

Manufactured in the UK

OptiGene Ltd.,
Unit 5 Blatchford Road, Horsham,
West Sussex, RH13 5QR, UK

Tel: +44 (0) 1403 274980

FAX: +44 (0) 1403 271017

Notice to purchaser: limited license

The purchase price of this product includes a limited, non-transferable license under European Patents EP1,020,534, EP1,873,260, EP2,045,337, EP1,416,055, EP1,724,361 or their foreign counterparts, owned by Eiken Chemical Co., Ltd.

Version 3.0 (May 2022)